


## Instructions for using Secure-Seal™ Imaging Spacers

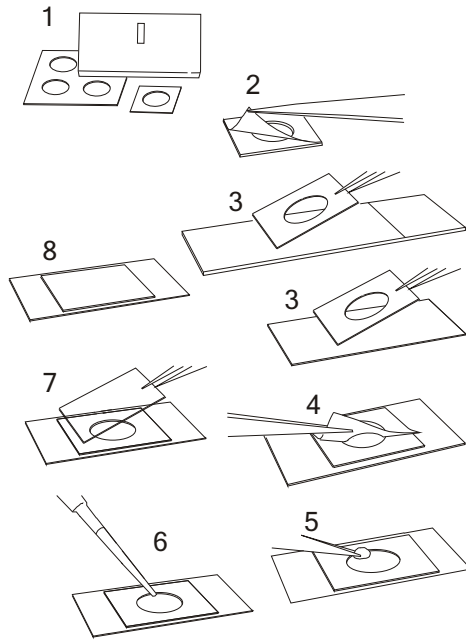

1. Trim to size using a razor blade or scissors (if desired).
2. Using forceps, remove adhesive liner from one surface.
3. Apply the spacer, adhesive side down onto the surface of a glass slide or coverslip. Press gently to seal.
4. Remove the other adhesive liner.  
*Note: To increase chamber depth add multiple spacers.*
5. Add specimen.
6. Add mounting medium.
7. Coverslip (if desired), applying gentle pressure over the adhesive area to seal.
8. Image