


## ProCrystal™ Protein Crystalization Covers


These instructions depict use of item 845236 which is pre-cut and has a paper release liner; however, they apply also to those pre-cut products that have a clear plastic liner.

1. Place a plate and cover onto a work surface with the cover adhesive facing up (Figure 1).

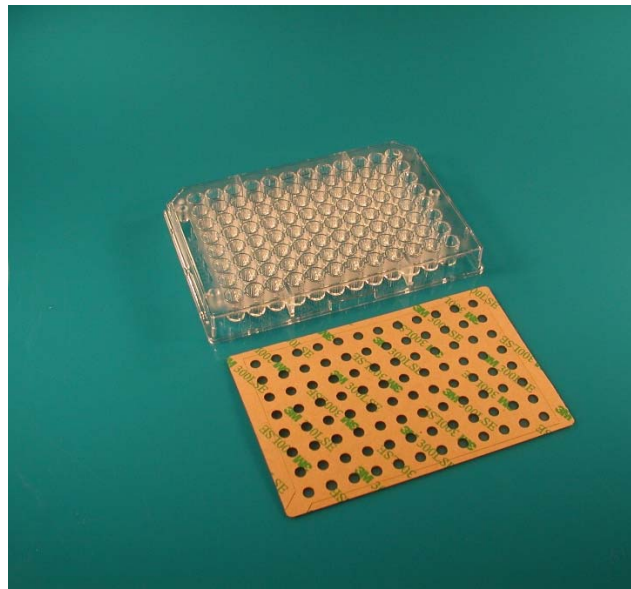


**Figure 1**

2. Expose the adhesive by peeling the paper release liner off of the cover (Figure 2).

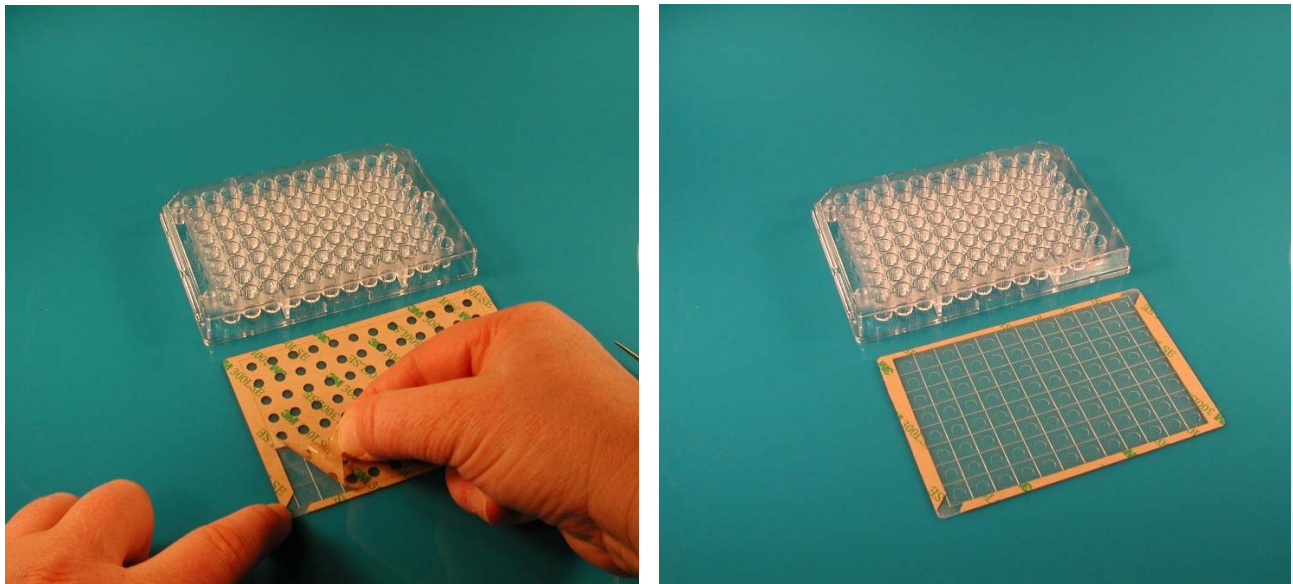


**Figure 2**

3. Holding the protected edges of the cover as shown in Figure 3, place it adhesive side down onto the plate, aligning the cover to the plate over the wells.

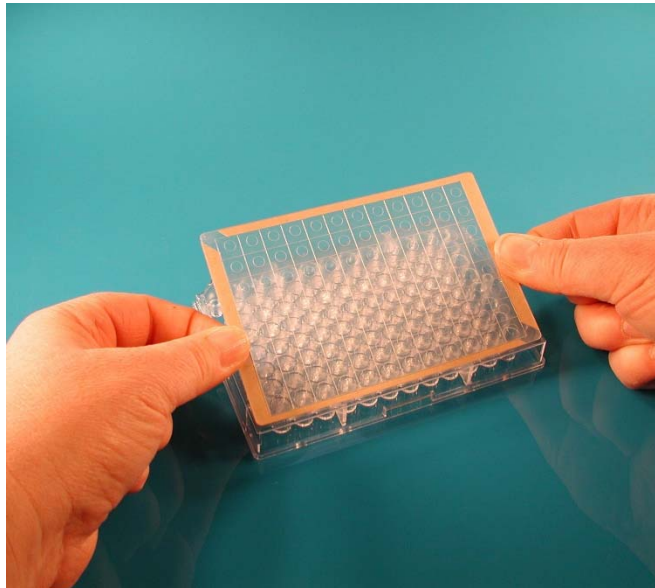


**Figure 3**

4. Press firmly to seal the cover to the wells as shown in Figure 4.

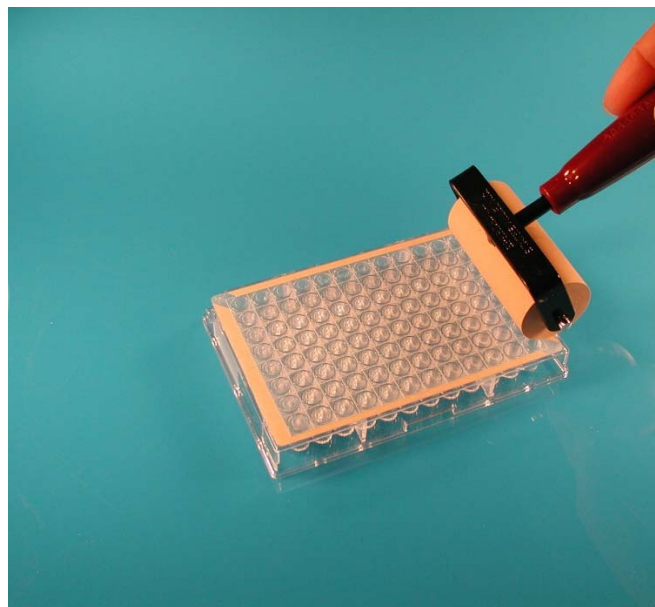


**Figure 4**

5. Remove the backing liner and frame by peeling it slowly away from the plate (Figure 5).

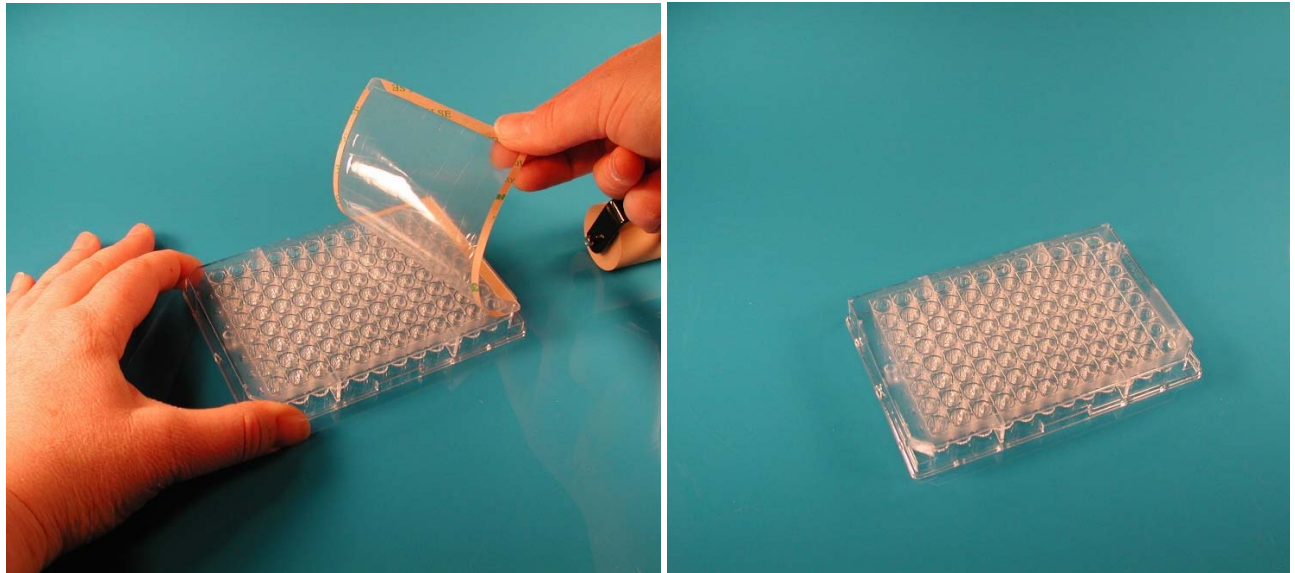

**Figure 5**


6. Remove an individual well cover by inserting a forceps into the corner of the well cover. Grasp the cover firmly and peel it off as shown in Figure 6.

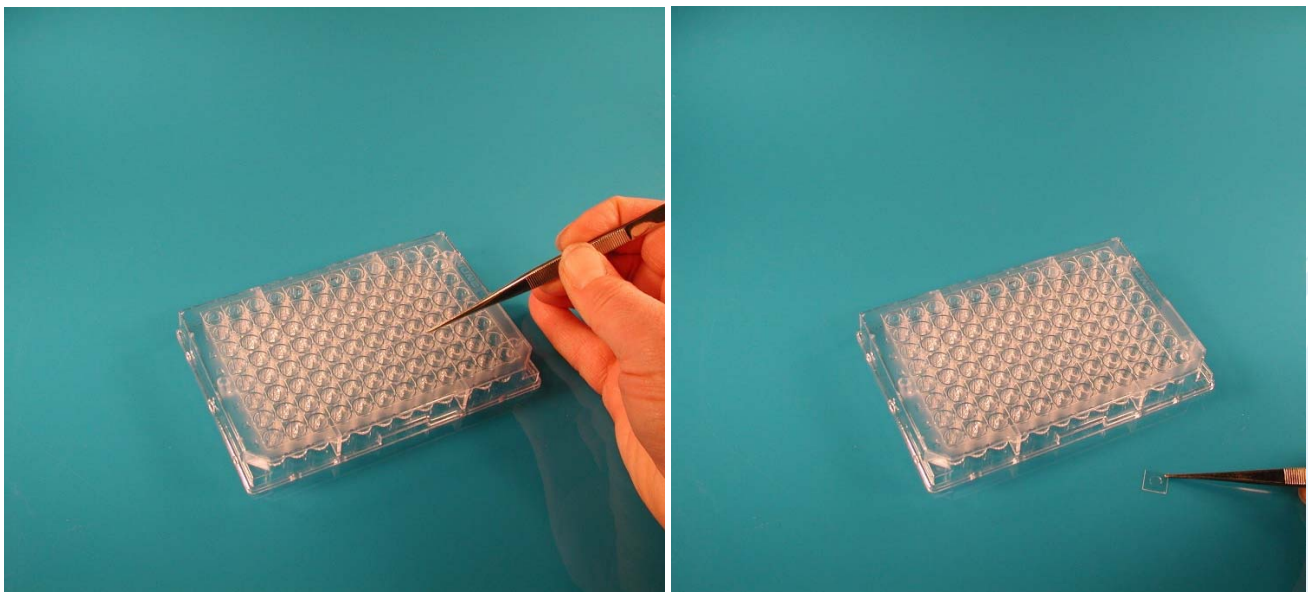

**Figure 6**

ProCrystal™ Protein Crystallization Covers

7. Figure 7 is a close up view of the well cover.

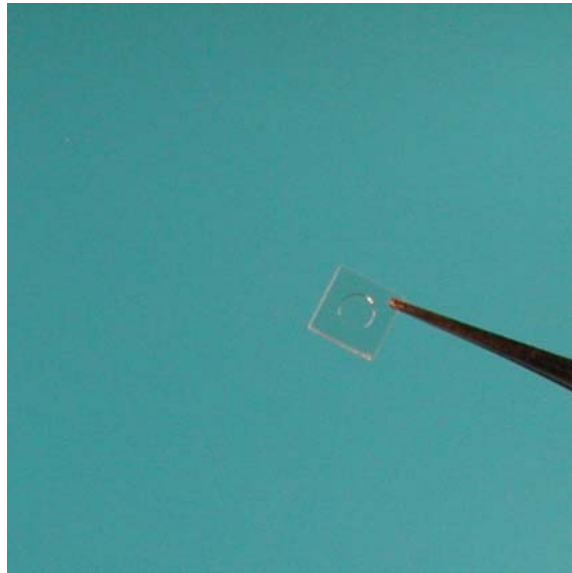

**Figure 7**