



## APPLICATION

1. Trim to size using a razor blade or scissors (if desired) as shown in figure 1.

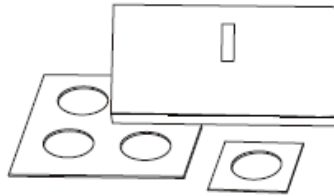


**Figure 1**

2. Use a forceps to remove the adhesive liner from one surface as shown in Figure 2.

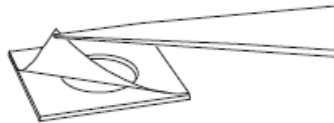

**Figure 2**


3. Apply the spacer, adhesive side down onto the surface of a glass slide or coverslip as shown in Figure 3. Press gently to seal.

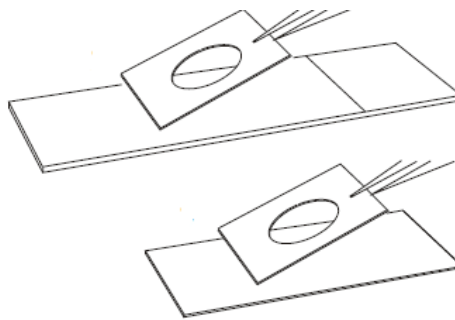

**Figure 3**


4. Remove the other adhesive liner as shown in Figure 4.

*Note: To increase chamber depth, add multiple spacers.*

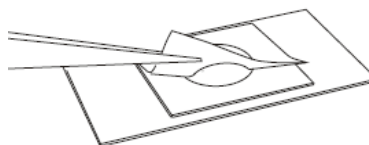


**Figure 4**

5. Add the specimen as shown in Figure 5.

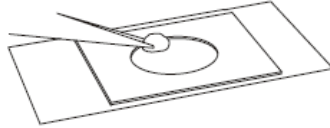


**Figure 5**

6. Add the mounting medium as shown in Figure 6.

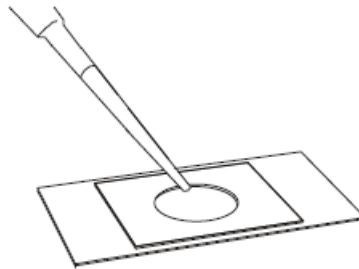

**Figure 6**

7. Add a coverslip (if desired) as shown in figure 7. Apply gentle pressure over the adhesive area to seal.

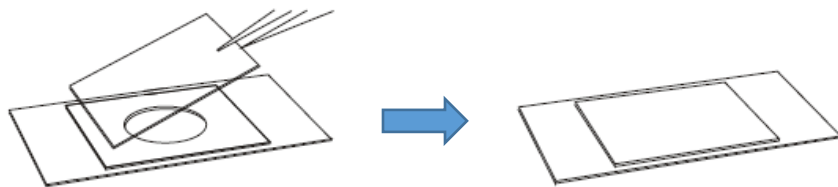

**Figure 7**

8. Image